


Jumbo Pipettes

General

The Jumbo Pipette is a single use, disposable, device for sampling low and medium viscosity liquids.


Jumbo Pipettes are assembled in a medical grade

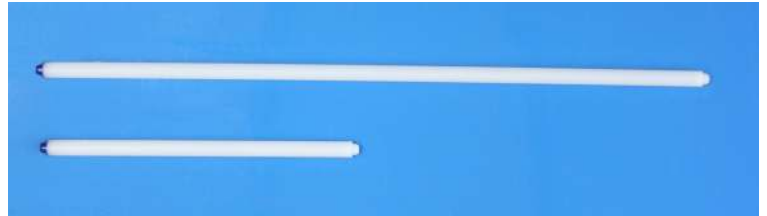

Operation

①

Hold the BLUE end.
Press your thumb onto tube until you are at the sampling depth

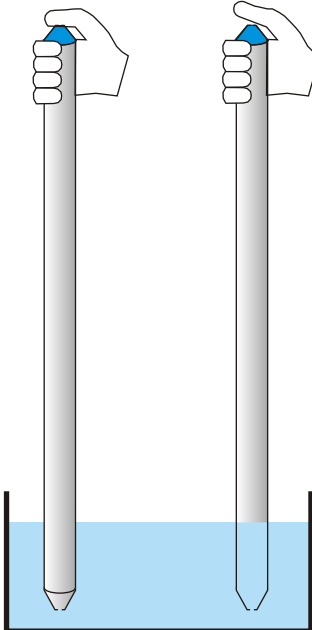


②

Raise your thumb until the levels equalise



③

Press thumb onto tube to seal in the sample

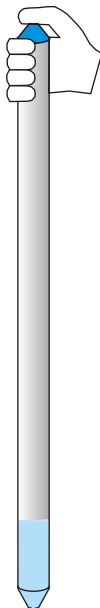

Cleaning

1. This is a single use device - do not clean
2. Do not re-use
3. Do not remove the top or bottom tapered fittings - they are welded in place and so can will not be secure when replaced

If you have any questions regarding the use of this device please contact Sampling Systems